

# Atom Physics

## Instadose App Setup



1. Open your preferred app store on your mobile device and search for the 'Instadose' app. Select the download button.
2. Your device's Bluetooth must be turned on to transfer data through the app.
3. Open the app and enter your username and password.  
**Note: If you forget your username or password, you can reset it or contact Atom Physics.**
4. A welcome screen will appear notifying you of any badge reports. Click the up arrow in the bottom right corner to continue to the app.



Shown above: the login screen, welcome screen, badges tab, and badge details

For help with the Instadose app, contact Atom Physics at 303.748.5499 or [info@atomphysics.com](mailto:info@atomphysics.com)

# Atom Physics

## Instadose Badge Reading



Your badge has already been initialized and is constantly collecting data. However, logging readings in the app is the only way to see the dose exposure readings collected by the badge. It is recommended to do this monthly.

1. Open the instadose app (with Bluetooth on) and select the 'Badge Reader' icon on the bottom of the screen
2. Hold the badge near your device. Press and hold the button on the back of the badge for 5 seconds until the indicator light stays green for one second.  
**Note: If a red light appears, wait one minute then try again. If the problem persists, contact Atom Physics at 303.748.5499 or [info@atomphysics.com](mailto:info@atomphysics.com) for troubleshooting.**
3. Wait up to 60 seconds for all of the dose readings to transfer to your device. The indicator light will stay green for one second when it is complete. Wait for the 'Finished Reading' screen to confirm the data transfer.

**Note: An asterisk (\*) denotes a reading lower than 1 mrem**

